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Prospekt výrobce

don whitley
scientific
excellence in microbiology

Whitley Workstations Whitley A35 Anaerobic Workstation

2

Sample manipulation within a secure anaerobic environment is now as easy as working aerobically at the laboratory bench. The Whitley A35 Anaerobic Workstation has been created by listening to customer needs and fulfils the diverse specifications of both research and clinical customers.

Even with a relatively small footprint incubation capacity is maximised and creates the perfect controlled environment, for large or small numbers of samples.

Complete with a built-in, rapid airlock and an intuitive touchscreen control panel, this workstation also allows you the option of working completely glove and sleeve free if fitted with instant access porthole system. You can, however, specify your A35 with a conventional sleeve and cuff porthole system.

Both the working and storage areas of the chamber are temperature controlled with an impressively low temperature gradient throughout. This allows the full internal chamber volume to be used for incubation if desired.

The optional Instant Access Porthole System eliminates gloves and sleeves - uses eight times less gas than conventional porthole systems

Rapid, built-in airlock with mechanical and electrical interlocks - no risk of compromising conditions inside the chamber, flushed with nitrogen as an additional cost saving

Innovative touch screen interface with password-protected user levels

Available with Removable Front

Dual gas operation

Automatic commissioning cycle

Automatic de-humidity control system – no user maintenance necessary

Available with integrated Anaerobic Indicator System

13

www.dwscientific.co.uk

Prospekt výrobce

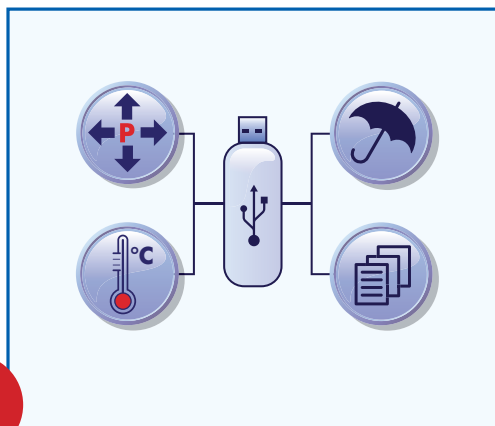
Easily removable front for equipment transfer and thorough cleaning

9cm Front Loading or 15cm Side Entry Letterbox - Ideal for introducing individual samples quickly.

Optional internal country-specific mains power sockets.

Convenient USB memory stick download of temperature, humidity, internal pressure and events log.

Intuitive operational software accessed via a full colour touch screen interface

The new directional workstation spotlight assists in specimen examination and the early identification of sample growth.

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Prospekt výrobce Whitley A35 Anaerobic Workstation fitted with the Instant Access Porthole System

The integral 12 litre, rapid-cycle airlock has the capacity to introduce up to 40 Petri dishes or small items of equipment into the workstation in 60 seconds.

An optional 9cm or 15cm letterbox can be specified, ideal for quickly transferring all sizes of dishes and small items.

Factory-Fitted Options

A07208 9cm Front Loading Letterbox - Ideal for quickly introducing small quantities of samples.

A06103 15cm Side Entry Letterbox - Ideal for quickly introducing small quantities of samples.

A06106 Double Internal Socket - (1 Amp maximum)

A02945 Spare Cable Gland - The standard cable gland allows cables and probes (3mm - 7mm in diameter) to be introduced into the chamber without compromising internal conditions. More than one gland can be fitted. Glands to accommodate larger diameters are available please discuss with our technical staff.

A07201 Anaerobic Indicator System (includes data logging) - Provides confirmation that anaerobic conditions exist inside the workstation and an early indication if conditions begin to vary.

14

9

Software/Accessories

A07200 Data Logging - For temperature, humidity, chamber pressure and events log.

A06110 Trolley - Custom-designed trolley frees up bench space and allows the workstation to be moved easily.

A06114 Internal Storage Tray / maximum of 2 per unit - One or two trays can be positioned under the shelf. This is either a factory-fitted option or trays can be fitted on-site.

A00006 Gas Leak Detector

A03872 10 Plate Petri Dish Rack Coloured Label Set (x140 / 7 different colours) - Enables users to colour code individual stacks of Petri dishes. Notes can be written on the reuseable tags using either water soluble or permanent markers.

A02534 10 Plate Petri Dish Rack

A07211 A35/A45 Spotlight Assembly - assists in specimen examination and the early identification of sample growth.

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Prospekt výrobce

A35 Instant Access Portholes - just three easy steps to use

About Don Whitley Scientific

For over thirty years we have pioneered the development and use of modified atmosphere workstations for microbiology and cell culture applications

More than 2000 of our workstations are in use in fifty countries

Worldwide technical support

Scientific guidance available from our own in-house contract laboratories

A major advantage of the sleeveless system is that it uses eight times less gas when entering the workstation compared with a conventional porthole system. The airlock is flushed with nitrogen as an additional cost-saving.

If you need to regularly deep clean your workstation chamber or want the facility to introduce small items of equipment or samples in bulk, you may prefer to order your A35 with a removable front. A removable front can be specified with Instant Access Ports or with a conventional port system using traditional sleeves and cuffs. The conventional port option is ideal for users wishing to enter the workstation less frequently for longer working periods.

The A35 is also now available with a fully integrated Anaerobic Indicator System. This factory-fitted option displays real time oxygen levels using coloured segments - green, amber or red - depending upon the percentage of oxygen present. The data gathered can be downloaded for further analysis.

Whatever your application, we can tailor the Whitley A35 Workstation to your requirements. For example, if you want to be able to use electrical equipment inside the chamber, you could include the option of an internal double electrical socket.

The A35 Workstation will accommodate between 400-600 x 90mm Petri dishes depending on whether plate carriers are used and which accessories and system options have been incorporated.

Specification

1

A07000	230V	Whitley A35 Anaerobic Workstation External dimensions (L x H x D) 1255mm x 710mm x 720mm (without removable front)
A07001	115V	

3

In the interests of a policy of continuous product improvement the company reserves the right to alter product specifications without prior notice. All rights reserved. © 2012 Don Whitley Scientific Limited.

Don Whitley Scientific Limited

14 Otley Road, Shipley, West Yorkshire, BD17 7SE, England.

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Prospekt výrobce

Installation Requirements

Whitley A35 Anaerobic Workstation User Manual / Installation and Commissioning

Physical Requirements

If bench mounted, the bench allocated must be flat, level and of sufficient size to support the base of the instrument fully.

Height:	710mm	(28")
Length:	1248mm	(49.25")
Depth:	711mm	(28")
Weight	100kg	(220lbs)

Do not site the instrument near draughts caused by doors, windows or air conditioning systems.

Environmental Requirements

Location:	Indoor use only
Altitude:	Up to 2000m
Degree of Protection:	IP20
Atmosphere:	Absence of corrosive gas. Minimum dust.
Operating Temperature:	5°C above ambient to 45°C
Relative Humidity:	4 The set temperature must be at least 5°C above ambient for effective control of the chamber temperature.
	Maximum 80% RH, without significant condensation, for temperatures up to 31°C, decreasing linearly to 50% RH at 40°C.
Applicable Rated Pollution Degree:	Pollution degree 2

Maintaining Good Anaerobic Conditions

Good anaerobic conditions are dependent upon several physical and environmental factors, which are summarised in the diagram below. When these factors combine and interact as intended, laboratory personnel can successfully and repeatedly cultivate anaerobic organisms.

If the conditions inside the workstation are not anaerobic, some bacteria will not grow well. This does not mean that anaerobic bacteria are not present.

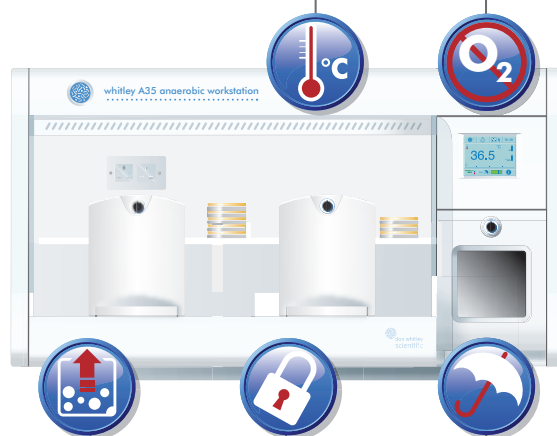
Stable Temperature

Temperature distribution within the workstation should be uniform. DWS workstations maintain temperature within $\pm 0.6^{\circ}\text{C}$ of the set point across the working and incubation areas of the chamber.

7

No Oxygen

Anaerobic gas mixture contains 10% hydrogen, 10% carbon dioxide and 80% nitrogen. Residual oxygen is removed by combining with hydrogen in the presence of a catalyst to produce water.

Catalyst + Anotox™

Catalyst and Anotox must be in good condition and not allowed to get wet. The typical life expectancy of catalyst is between 6-12 months depending on the types of bacteria being grown. Anotox is essential for maintaining ideal internal conditions by absorbing volatile fatty acids and hydrogen sulphide, both of which can be byproducts of the growth of anaerobic bacteria. Anotox has been shown to substantially prolong catalyst life. Don Whitley Scientific recommends that catalyst and Anotox are replaced together.

10

Relative Humidity Control

To maintain ideal bacterial growth conditions, agar must not be allowed to dry out during sample incubation. Excessive internal moisture may promote bacterial cross contamination.

A Sealed Workstation Chamber

Workstations are operated at a slight positive pressure so that any small leaks allow anaerobic gas to leak out, instead of allowing oxygen to enter the workstation.

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Whitley A35 Workstation

Enter Keyword
whitley-
scientific-
limited)

OVERVIEW SPECIFICATIONS ACCESSORIES IMAGES/VIDEOS DOWNLOADS TESTIMONIALS GOT A QUESTION? REQUEST A CALLBACK

Specifications


Capacity (Litres)	300
Capacity (Petri Dishes)	400 / 600
Port/Airlock Capacity	40 plates via airlock
Porthole System	Manual or Instant Access Ports
Gas Supplies	ANO2/N2
Footswitch	Wireless
Auto Sleeve Gassing	Option available
Internal Mains Socket	Option available
Spotlight / Inspection Lamp	Option available
Single Sample Entry	Option available
ANO2 Conditions Monitor	Option available
Refrigeration	Not applicable
Data Logging	Option available
Airlock Cycle Time (seconds)	60
Extra Cable Glands	Option available
Microscope	Contact us for further information
Automatic Dehumidifier	Fitted as standard
Automatic Humidifier	Not applicable
Removable Front	Option Available
Trolley	Option Available
Dimensions (w/d/h - mm)	1255 / 720 / 710
Weight (lbs/kg)	220 / 100
Temperature Range (°C)	5°C above ambient up to 45°C
Touchscreen control	Yes

15 cm Single Plate Entry System on the A35

Rollover Image to Zoom

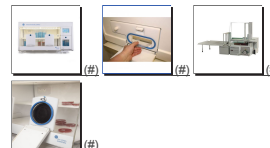ADDITIONAL IMAGES & VIDEOS

View further images & videos (#)

PRODUCTS

Whitley Anaerobic Workstations
(<http://www.dwsscientific.co.uk/whitley-anaerobic-workstations/>)

Whitley Hypoxic Workstations
(<http://www.dwsscientific.co.uk/whitley-hypoxic-workstations/>)

Whitley HEPA Workstations
(<http://www.dwsscientific.co.uk/whitley-hepa-workstations/>)

Whitley Instrument Workstations
(<http://www.dwsscientific.co.uk/whitley-instrument-workstations/>)

Whitley Microaerophilic Workstations
(<http://www.dwsscientific.co.uk/whitley-microaerophilic-workstations/>)

Spiral Plating
(<http://www.dwsscientific.co.uk/spiral-plating/>)

Lab Automation
(<http://www.dwsscientific.co.uk/lab-automation/>)

Tissue Culture Systems
(<http://www.dwsscientific.co.uk/tissue-culture-systems/>)

More Products
(<http://www.dwsscientific.co.uk/more-products/>)

ABOUT US

About Us
(<http://www.dwsscientific.co.uk/about-us/>)

Certificates
(<http://www.dwsscientific.co.uk/certificates/>)

Testimonials
(<http://www.dwsscientific.co.uk/testimonials/>)

Export Sales
(<http://www.dwsscientific.co.uk/export-sales/>)

History
(<http://www.dwsscientific.co.uk/history/>)

Products & Services
(<http://www.dwsscientific.co.uk/products-and-services/>)

Overseas Distributors
(<http://www.dwsscientific.co.uk/overseas-distributors/>)

Exhibitions
(<http://www.dwsscientific.co.uk/exhibitions/>)

Equipment Rental Example
(<http://www.dwsscientific.co.uk/equipment-rental-example/>)

Vacancies Team
(<http://www.dwsscientific.co.uk/vacancies-team/>)

Workstation Poster Grant
(<http://www.dwsscientific.co.uk/workstation-poster-grant/>)

SERVICE & ACCREDITATION

Service & Accreditation
(<http://www.dwsscientific.co.uk/service-and-accreditation/>)

UKAS Calibration / Validation
(<http://www.dwsscientific.co.uk/ukas-calibration-validation/>)

Contract Laboratory
(<http://www.dwsscientific.co.uk/contract-laboratory/>)

RECEIVE OUR EMAIL UPDATES

To receive our latest news, products & exhibitions

enter email address

SUBMIT

© 2015 Don Whitley Scientific Limited
Registered Office: 14 Otley Road, Shipley, West Yorkshire, BD17 7SE, UK. Registered in England No. 1342672
Terms and Conditions (<http://www.dwsscientific.co.uk/terms/>) | Privacy Policy (<http://www.dwsscientific.co.uk/privacy/>) |
Cookie Policy (<http://www.dwsscientific.co.uk/cookies/>) | Site Map (<http://www.dwsscientific.co.uk/sitemap/>)
(<https://plus.google.com/117563081616039026978>)

Web Design (<http://www.dwsscientific.co.uk/web-design/>) - Rejuvenate Productions

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Prospekt výrobce

The A35 Hypoxystation

Whitley A35 Anaerobic Workstation User Manual / Introduction

The A35 Anaerobic Workstation

WHITLEY A35 ANAEROBIC WORKSTATION

MANUFACTURED BY
DON WHITLEY SCIENTIFIC LIMITED
WEST YORKSHIRE, BD17 7SE, UNITED KINGDOM

T: +44 (0)1274 595728 F: +44 (0)1274 531197
www.dwscientific.co.uk

WARNING

INSTALLATION AND COMMISSIONING SHOULD ONLY BE UNDERTAKEN BY PERSONNEL WHO HAVE RECEIVED PRODUCT TRAINING FROM DON WHITLEY SCIENTIFIC LIMITED

INPUT PRESSURES: Minimum 4 bar (60 psi) / Maximum 6 bar (90 psi)

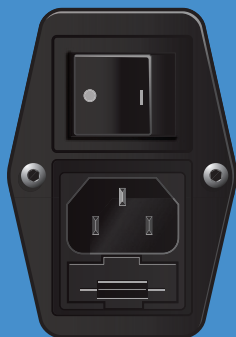
NITROGEN

ANAEROBIC GAS

Hydrogen 10%
Carbon Dioxide 10%
Nitrogen 80%

POWER ENTRY MODULE

DISCONNECT MAINS POWER SUPPLY BEFORE REPLACING FUSES F1 AND F2

F1 T10A
F2 T10A

Rear Panel

Positioned on the rear of the instrument behind the airlock.

This panel contains gas and electrical connections, fuses, manufacturers contact details and other essential information.